

Liquid for inhibiting algae growth

Supplied in a 1 liter dosing bottle, **Clinet[®] anti-algae** is used, once diluted to 0.1%, to prevent algae growth in laboratory water baths.

Application

Algae-inhibiting liquid for laboratory water baths.


How to use

Dilute **Clinet[®] anti-algae** to 0.1%. Pour a 20 ml dose into a water bath of 20 liters of distilled water.

Renew the dose every week. Change the water every month.

Presentation

1 liter bottle with dosing dispenser.

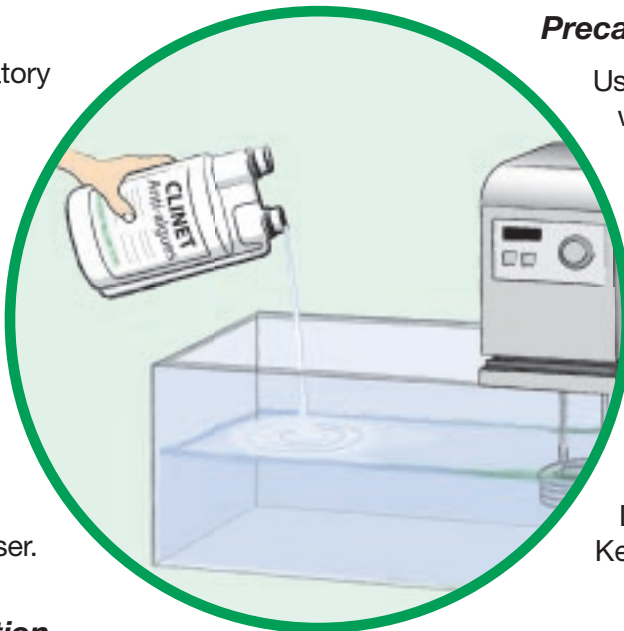

Precautions

Use only demineralized or deionized water, as the algicide activity is reduced by hard water.


Never mix with anion-active detergents and/or strong oxydants (such as bleach...).

In case of contact with the eyes, rinse abundantly with water and consult a doctor.

Do not swallow.
Keep out of reach of children.

Composition

Quaternary ammonium, forms a complex with calcium, preservative.

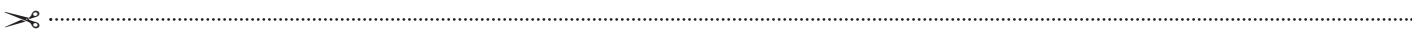

- Please send me :
- a price offer for Clinet[®] anti-algae.
 - technical information on Clinet[®] anti-algae.
 - a free sample of Clinet[®] anti-algae.

Company

Name

Position

Address

Zip code/City

Country

Tel. Fax

E-mail

Hygiene & Security

	information	samples
Clinet[®] 101 (chlorhexidine cleaning wipes)	<input type="checkbox"/>	<input type="checkbox"/>
Clinet[®] 202 (indivi. chlorhexidine cleaning wipes)	<input type="checkbox"/>	<input type="checkbox"/>
Clinet[®] 303 (bactericidal spray)	<input type="checkbox"/>	<input type="checkbox"/>
Anabac[®] (autoclave deodorant)	<input type="checkbox"/>	<input type="checkbox"/>
BioParfum[®] (odor neutralizers)	<input type="checkbox"/>	<input type="checkbox"/>

BagSystem[®] for sample preparation

BagMixer[®] (sample blenders)	<input type="checkbox"/>	
BagFilter[®] (sterile blender bags)	<input type="checkbox"/>	<input type="checkbox"/>
BagPipet[®] (automatic pipette)	<input type="checkbox"/>	

Please fax this request to your local dealer or to :

www.ert.pt
info@ert.pt

Tel. 217 142 066 - Fax. 217 145 674

Rua Soeiro Pereira Gomes, Lote 5, Cave Dt.
2845-387 Amora Portugal